



## Pocket Acidity Meter

# PAL-ACID2

Cat.No.4652

The PAL-ACID2 measures the total acidity in sample and converts it into tartaric acid concentration.

### How to Measure

Pipette 0.1ml (100 $\mu$ l) of the sample and dispense it into the quinone reagent solution.

Fill the sample stage with the solution, put the metal anti-volatile cover on, and then press the START key.

The acid content is displayed in grams per liter.

### Pocket Acidity Meter PAL-ACID2 Test Kit



The kit comes complete with 10 bottles of quinone reagent solutions (5ml each), a micropipette, 10 micropipette tips, 10 disposable plastic pipettes, 1 bottle of calibration solution, an anti-volatile cover, and 2 AAA batteries.

### Consumable Supplies

- Micropipette **RE-79401**
- Calibration solution **RE-130002**
- Quinone reagent solution **RE-99432** 10 pcs
- Quinone reagent solution **RE-99430** 20 pcs
- Quinone reagent solution **RE-99431** 50 pcs

One pipette tip per bottle is included.

### Commonly Measured Samples

	(g/l)
Grapes (Pione)	6.0
Grapes (Delaware)	5.9
Grapes (Kyoho)	4.7
Red wine (Sangiovese)	5.3
Red wine (Cabernet Sauvignon)	4.9~6.5
White wine (Chardonnay)	5.0~7.3
Grape juice	2.8~5.4

PAL-ACID2 Test Kit		Cat.No.4652
Measurement Range	Tartaric acid concentration (total-Tartaric acidity conversion [g/l]) 1.0 to 40.0 (g/l)	
Repeatability	1.0 to 20.0 (g/l) $\pm 0.5$ (g/l) 20.1 to 40.0 (g/l) $\pm 1.0$ (g/l)	
Measurement Temperature	10 to 40°C (Automatic Temperature Compensation)	
Standard Accessories	<ul style="list-style-type: none"> <li>• Micropipette : RE-79401</li> <li>• Calibration solution : RE-130002</li> <li>• Quinone reagent solution 10 pcs :RE-99432</li> </ul>	
Power Supply	2 x AAA batteries	
Minimum Indication	0.1 (g/l)	
Ambient Temperature	10 to 40°C	
Dimensions & Weight	55(W) $\times$ 31(D) $\times$ 109(H)mm, 100g (Main Unit only)	



**Video manual available!**

For more details, visit us at [www.atago.net/manual](http://www.atago.net/manual)

All ATAGO refractometers are designed and manufactured in Japan.

**HACCP GMP GLP**

ATAGO products comply with HACCP, GMP, and GLP standards.